



Continuous Dynamic Analysis and Quasi-Static Measurement of Spider Silks

Application Note

John Swindeman
Agilent Technologies



Figure 1. Forcible silking of *Argiope argentata*.

Introduction

Nanoscale characterization of biological materials continues to be a challenge in the fields of mechanics and material testing. One example of this type of research is the characterization of the mechanical properties of spider silks. Such silks have an astounding strength-to-weight ratio that has attracted the interest of a number of communities, ranging from medical to military.

Complexities associated with spider silk research include the small and hard-to-measure diameter of the silks, difficulties harvesting and gripping the samples, and the ability to obtain meaningful quasi-static results over large strain ranges as well as to measure various dynamic characteristics.

Sample Preparation

Among the greatest challenges of measuring such small fibers are harvesting the specimens and determining a suitable method for mounting them. The samples can be

harvested in several ways. Silks can be collected from native sources such as webs, egg sacs, draglines, and the swathing silks used to wrap prey. As shown in Figure 1, silks can also be harvested via forcible silking, a process that entails pulling fibers directly from the spinnerets of anesthetized spiders.

Scientists at Agilent Technologies used a template technique to prepare spider silk samples of similar size. A circular or rectangular window was cut into the template and a specimen was glued across the window. The template was then clamped at both ends in an Agilent T150 universal testing machine (UTM). Next, the template was cut on either end of the window so the material could be tested.

Displacement Sensing

A number of spider silk samples were tested using the T150 UTM. The three spider silks presented in this application note are from the species *Uloborus diversus*, *Argiope argentata* (silver garden spider), and *Latrodectus*

hesperus (western black widow). Figure 2 and Figure 3 show an SEM image and quasi-static results for *Uloborus diversus*. This sample is of particular interest, as it is only 500 nm in diameter. These fibers can only be seen with the naked eye by the reflection of strong light. Note that the load portion of the curve in Figure 3 is in micronewtons.

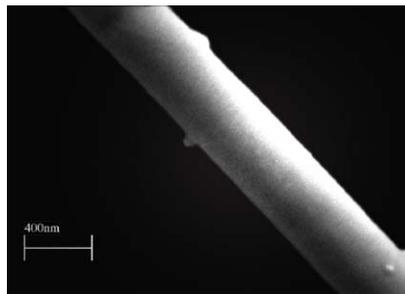


Figure 2. SEM image of 500 nm *Uloborus diversus* spider silk.

Uloborus diversus has some notable material characteristics, including an ultimate tensile strength of 1.150 GPa and a modulus as high as 13.861 GPa. Furthermore, a sample with a stiffness as low as that shown in Figure 3 (~0.124 N/m) presents significant challenges and exemplifies the capabilities of the T150 UTM.

In Figure 4, the engineering stress/strain curve can be seen for silk harvested from the species *Argiope argentata*. This silk is often referred to as a “sticky silk” and is not as dry as the other two silks mentioned above. An interesting characteristic of this silk that differs from the other two is the extreme strain that it can withstand. The stress/strain curve is very shallow out to nearly 300% strain before it starts to rise. In fact, it was pulled to nearly 430% strain before it broke. This remarkable behavior is what allows this material to perform its function so well.

Other material properties of interest include the dynamic characteristics (storage and loss modulus) and how these properties evolve as the static strain in the sample is increased. This is done by continuously superimposing an oscillatory force on the loading of the specimen. From this oscillating force and the resulting displacement oscillation, meaningful dynamic data can be obtained. The ability to make such measurements as the static strain progresses to very large values is a unique capability of the T150 UTM.

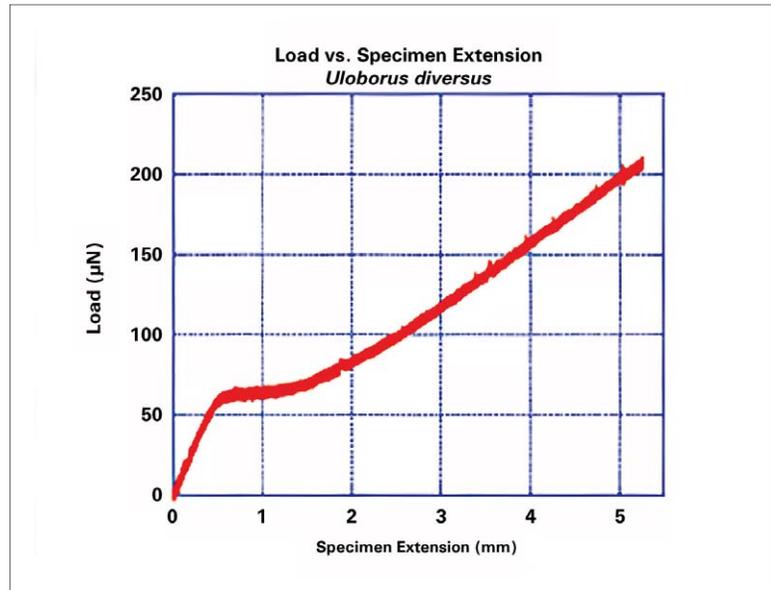


Figure 3. Loading curve for *Uloborus diversus*.

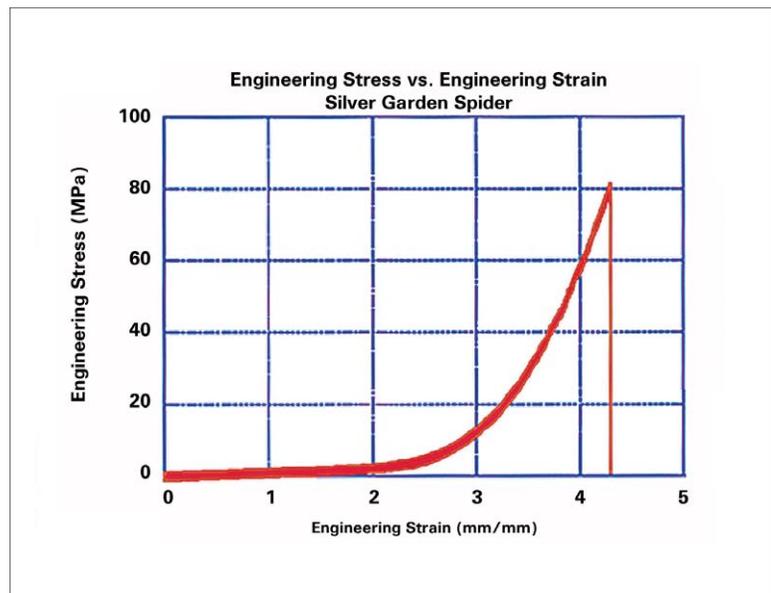


Figure 4. Loading curve for *Argiope argentata*.

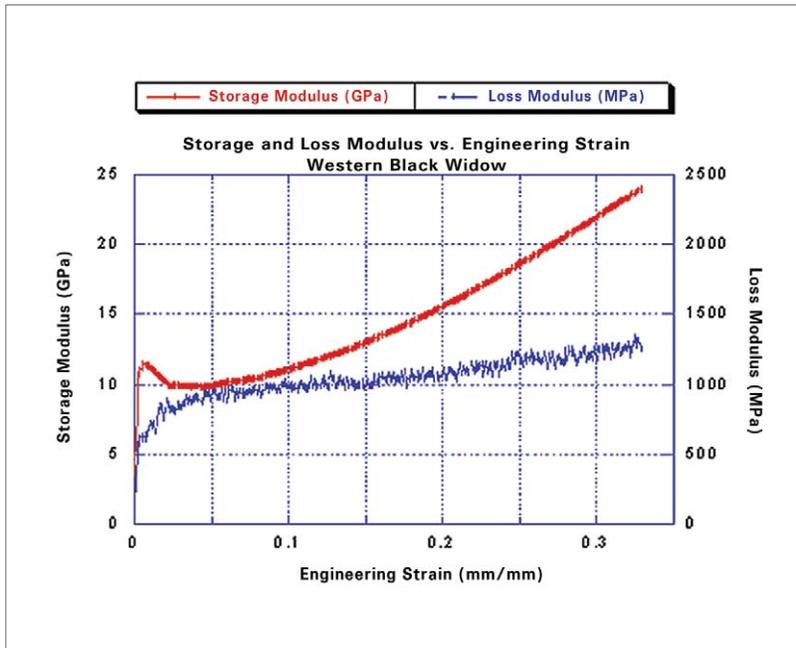


Figure 5. Storage and loss modulus for *Latrodectus hesperus* (western black widow).

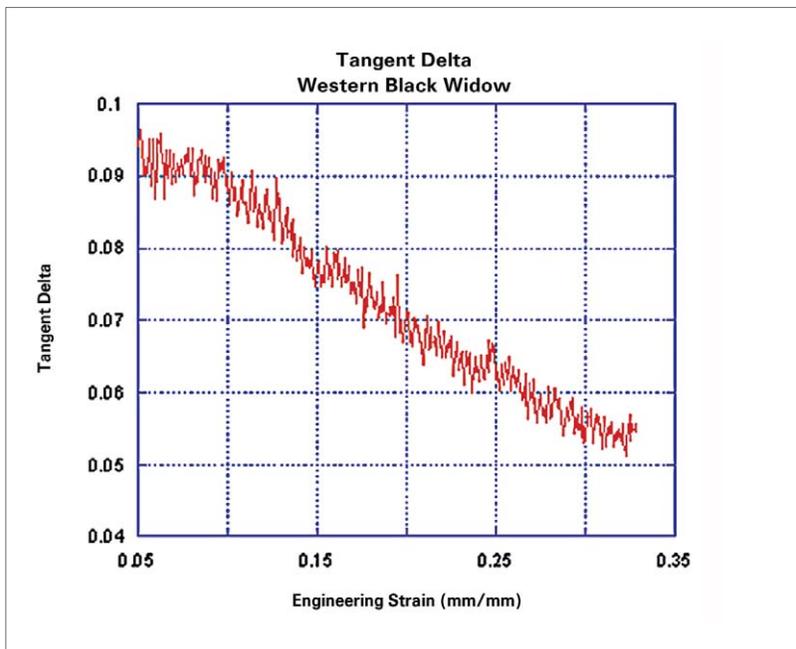


Figure 6. Tangent delta for *Latrodectus hesperus*.

Figure 5 and Figure 6 show the storage modulus, loss modulus, and tangent delta respectively for the spider species *Latrodectus hesperus*. Tangent delta is the ratio of loss modulus to storage modulus. The dynamic data is impressively quiet for a viscoelastic spider silk with a diameter of 5.574 microns. Accurate measurement of such properties offers valuable support for emerging research on spider silks.

Technology and Applications

The T150 UTM is a tensile testing machine that is capable of measuring force and extension on the nanoscale with exceptional noise control. Its extension axis, driven by a screw drivetrain, operates with better than 35 nm resolution using a rotary encoder; a nanomechanical actuating transducer head measures the force on the specimen with 50 nN resolution.

To enable mechanical properties to be determined continuously as the specimen is strained, the Agilent Continuous Dynamic Analysis (CDA) option allows the direct, accurate measurement of the specimen's stiffness at each point in the experiment. CDA makes it possible to determine storage and loss modulus, as well as to measure complex moduli over a range of frequencies.

Applications of the T150 UTM include yield of compliant fibers and biomaterials, dynamic studies of fibers and biomaterials, and tensile and compression studies of polymers.

Acknowledgments

Agilent Technologies would like to thank Todd Blackledge and Cheryl Hayashi of the University of California, Riverside for their insight.

Nano Mechanical Systems from Agilent Technologies

Agilent Technologies, the premier measurement company, offers high-precision, modular nano-measurement solutions for research, industry, and education. Exceptional worldwide support is provided by experienced application scientists and technical service personnel. Agilent's leading-edge R&D laboratories ensure the continued, timely introduction and optimization of innovative, easy-to-use nanomechanical system technologies.

www.agilent.com/find/nanoindenter

www.agilent.com

For more information on Agilent Technologies' products, applications or services, please contact your local Agilent office. The complete list is available at:

www.agilent.com/find/contactus

Phone or Fax

United States: (tel) 800 829 4444
(fax) 800 829 4433

Canada: (tel) 877 894 4414
(fax) 800 746 4866

China: (tel) 800 810 0189
(fax) 800 820 2816

Europe: (tel) 31 20 547 2111

Japan: (tel) (81) 426 56 7832
(fax) (81) 426 56 7840

Korea: (tel) (080) 769 0800
(fax) (080) 769 0900

Latin America: (tel) (305) 269 7500

Taiwan: (tel) 0800 047 866
(fax) 0800 286 331

Other Asia Pacific Countries:

tm_ap@agilent.com
(tel) (65) 6375 8100
(fax) (65) 6755 0042

Product specifications and descriptions in this document subject to change without notice.

© Agilent Technologies, Inc. 2009
Printed in USA, July 6, 2009
5990-4325EN



Agilent Technologies